

## va-Q-genic Handling Instruction

For va-Q-genic 2

**Note:** This product shall be stored and transported in upright position!

### Step I

(1) If sealing tape is applied, remove it or cut it open.  
Please use a safety cutter.

(2) Remove the tuck-in flap of the outer carton.



### Step II

Open the outer carton to reach the va-Q-genic 2.



### Step III

Open the top flap.

The outer package includes the va-Q-genic 2 and, if ordered, a data logger.

To prevent any damage to the dry shipper, please do not remove the dry shipper from its outer package.

**Note:** Further steps require appropriate personal protective equipment!



## Step IV

Remove the plug and the data logger, if included.



## Step V

Take out the two canisters one after the other.

Start with the centric canister because it keeps the eccentric canister in its position.

For detail see picture below, the difference between the two canisters is visible there.



## Step VI

Now the product can be placed inside the canisters.

The centric canister (1) and eccentric canister (2) are shown in the right picture.



## Step VII

Then proceed with the steps backwards to close the va-Q-genic 2.

Place the eccentric canister into its eccentric position, then the centric canister in the va-Q-genic 2.



## Step VIII

Put the plug back on the va-Q-genic 2 and place the data logger.

The arrow in the picture shows where to place the temperature probe.



## Step IX

---

Close the top flap.

If required, the va-Q-genic 2 can now be sealed or taped.



## Step X

---

Please follow the instructions on the outer carton to close it properly.

